



Octet Data Analysis HT Software Version 12

Newest features for advanced Epitope Binning and Quantitation analyses

Key features

- **NEW** BinChart for convenient binning of antibodies
- Advanced settings for fitting quantitation datasets
- Optimal visualization via overlaid fitted traces on top of raw quantitation binding traces
- **NEW** capability for reference subtraction within and across plates
- Optimized *Report Tab* for direct creation of exportable reports

Overview

The Octet® Data Analysis HT software is designed for fast analysis of large datasets from quantitation, kinetic and epitope binning assays. In version 12, the Data Analysis HT software offers faster and more advanced epitope binning and quantitation analysis options.

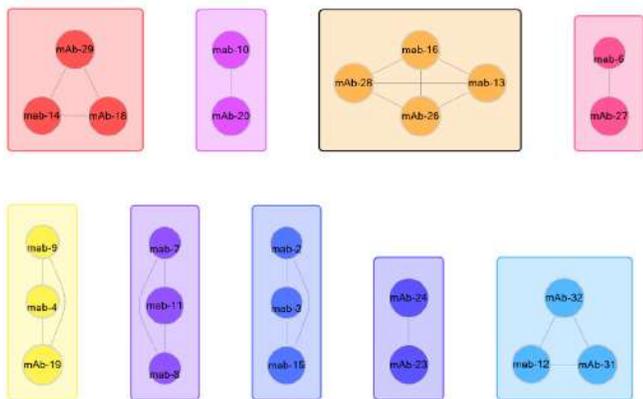


Figure 1: Example BinChart from a 32x64 screen. Each bin represents antibodies that block each other's ability to bind the antigen, and have similar blocking profiles when paired with other antibodies in the panel. This example was created using a low binding nm threshold setting, which resulted in fewer antibodies within each bin. When the mouse arrow is hovered over a bin, the top of the BinChart window also displays the list of antibodies within that bin.

BinChart

Epitope binning groups a panel of monoclonal antibodies (mAbs) into bins based upon similarities in antigen recognition sites or epitopes. This grouping is performed using cross competition assays, in which the competitive binding of antibody pairs to a specific antigen is characterized. If the binding of antigen to one mAb prevents the binding of another mAb, then these mAbs are considered to bind to similar or overlapping epitopes. If binding of a mAb to the antigen does not interfere with the binding of the other mAb, then they are considered to bind to distinct, non-overlapping epitopes. Two criteria must be fulfilled in order to assign mAbs into the same bin. First, all mAbs in the same bin should block each other's ability to bind the antigen. Second, all mAbs in the same bin should have similar blocking profiles when paired with other mAbs in the panel.

Version 12 introduces a BinChart feature which assigns screened antibodies to bins based on the binding threshold set during analysis. This binding threshold can be adjusted to analyze varying binning of antibodies as a result of the changing threshold (Figures 1 and 2).

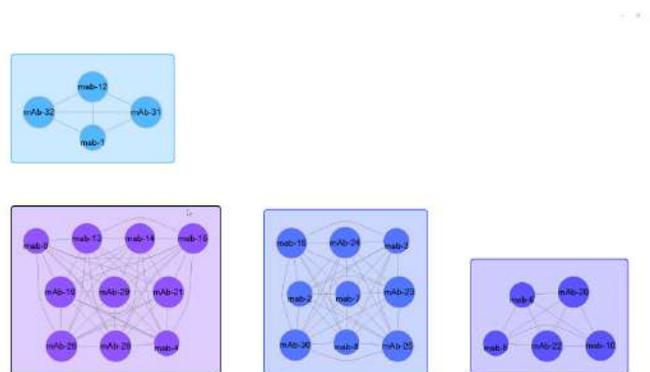


Figure 2: Example BinChart from a 32x64 screen. Each bin represents antibodies that block each other's ability to bind the antigen, and have similar blocking profiles when paired with other antibodies in the panel. In this example, a higher binding nm threshold setting was used, which resulted in more antibodies grouped in fewer number of bins compared to Fig 1.

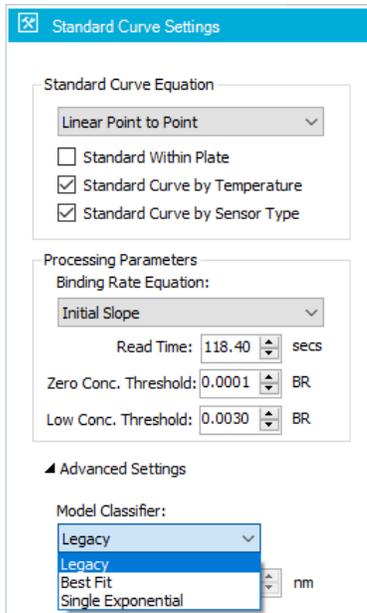


Figure 3: A screenshot of the Advanced Settings options for choosing among model classifiers.

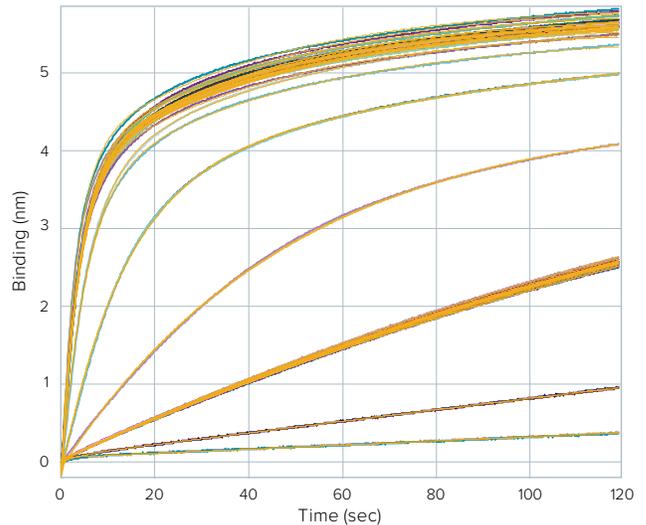


Figure 4: In Octet Data Analysis HT software version 12, fitted traces can now be overlaid on raw binding traces for optimal visualization. The *Best Fit* model classifier was used in this example.

Advanced quantitation

Version 12 adds new fitting model classifiers for advanced users. The model classifiers (Figure 3) help choose the best fit for the binding interactions. Version 12 also provides optimal visualization (Figure 4) by overlaying fitted traces on top of the raw binding traces.

- The *Legacy* model classifier chooses a fitting model (single exponential, double exponential) based on maximum slope of the raw binding traces.
- The *Best Fit* model classifier looks to minimize residuals – the difference between raw binding traces and fitted traces. If the difference is zero between different models (single exponential, double exponential), the *Best Fit* model classifier will choose the simpler of the two models.
- The *Single Exponential* model classifier forces fitting of raw binding traces with the single exponential mathematical model. This classifier might be best suited for known 1:1 binding interactions.

Report Tab

Take full advantage of the newly optimized Report Tab to create exportable reports with any number of data points. New features add capability to modify and sort data tables, and flexibility for font settings.

Ordering information

Part no.	Description
50-5029	SOFTWARE UPGRADE/OCTET SOFTWARE Upgrade of Octet instrument software to the latest version of software. Includes 4 seat licenses.
50-5043	SITE LICENSE UPGRADE - LATEST VERSION OF OCTET DATA ANALYSIS SOFTWARE. Upgrade of site license to the latest Octet standard software version at a physical site (single city or ZIP code) as defined in the purchase order. Requires previous purchase of Site License Octet Data Analysis Software.
50-5093	OCTET CFR SOFTWARE V12.X MODULE. Includes 1 seat license.
50-0296	OCTET CFR SOFTWARE V12.X VALIDATION PACKAGE



ForteBio
47661 Fremont Boulevard
Fremont, CA 94538
888.OCTET-75 or 650.322.1360
fortebio.info@moldev.com

ForteBio Analytics (Shanghai) Co., Ltd.
No. 88 Shang Ke Road
Zhangjiang Hi-tech Park
Shanghai, China 201210
salesops.china@moldev.com

Molecular Devices (UK) Ltd.
660-665 Eskdale
Winnersh Triangle
Wokingham, Berkshire
RG41 5TS, United Kingdom
+44 118 944 8000
uk@moldev.com

Molecular Devices (Germany) GmbH
Sauerbruchstr. 50
81377 München
Germany
+00800 665 32860