


Sartobind® Membrane Adsorbers, 4 and 8 mm Bed Height



### A Rising Star in Bioprocessing

The macroporous membrane structure and optimized design of the Sartobind membrane adsorbers enable efficient and ultra-rapid chromatographic purification. As a result, Sartobind is an economical solution for bioprocessing challenges and an asset for your purification platform.

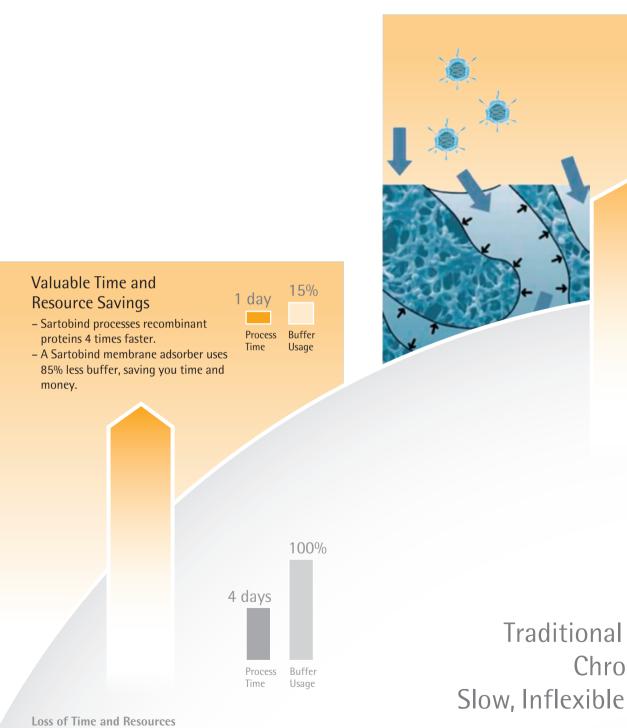

- Traditional column chromatography systems are slow in producing recombinant proteins.
- They waste considerable buffer resources.

Chro

# Discover the Favorable Performance and Economics of Sartobind Membrane Chromatography

#### **High Binding Capacity**

Sartobind membranes have a pore size of  $> 3 \mu m$ . This allows large proteins, bioparticles and viruses to enter the macro-porous membrane structure.

Consequently, Sartobind membrane adsorbers have a 10-times higher binding capacity for viruses, and about 200 times more polishing capacity than traditional chromatography columns.

#### Low Binding Capacity

Conventional column beads have a pore size < 100 nm. Small pores limit the access of large molecules to internal binding sites, resulting in a low binding capacity for large biomolecules, blood factors and viruses or virus-like particles. In addition, columns are operated at 10 to 30 times lower flow rates than are membrane adsorbers.

# Column matography: and Costly

# Compact Capsule Footprint Compared with a column or an ultracentrifuge, a Sartobind capsule is much more compact and easier to set up and use. As a

result, it requires less lab space and can be

conveniently and accurately scaled up.

 Disposable or reusable

No hardware investments



Large and Expensive Equipment



Pic. 1: Chromatographic columns have a limited loading capacity for polishing and virus capture. Furthermore, they are large and have a high investment cost.




Pic. 2: Likewise, expensive ultracentrifuges are difficult to scale and complex to operate.

# One Device for Capture and Polishing

The 8 mm capsules are void-volume-optimized, giving a sharp peak upon elution with minimal dilution. Compared to non-optimized capsules, this construction increases the dynamic binding capacity by at least one-third, and reduces the need for chromatography buffers by more than two-thirds.

#### **Optimized Design**

With conventional adsorbers, the extensive void volume has to be filled with buffer. Sartobind 8 mm capsules have a void-volume-minimizing core, creating an optimized flow channel, which ensures sharp elution peaks in bind and elute mode and less buffer consumption in the flow-through mode. The direction of flow is from the external upstream channel to the internal downstream channel.

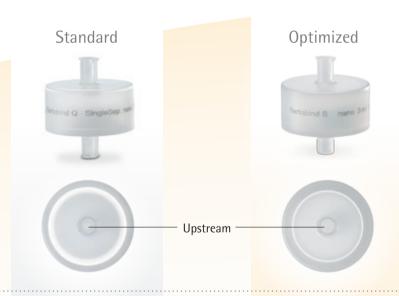

#### **Top View**

of the 3 mL version (with optimization of the flow channels)

## Cutaway of a 5" 70 mL Capsule


**Top View** 

zation)

of the 1 mL nano

version (no optimi-

Conventional adsorbers intended for polishing have no internal core. During operation, the open space is filled with buffer.

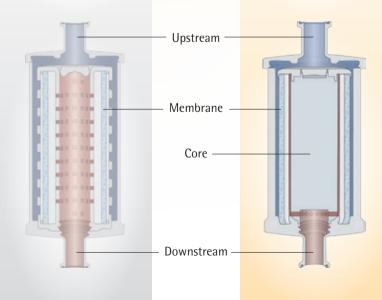


#### Cutaway of a 75 mL Capsule

The 8 mm capsule incorporates a core, which forms a small internal channel so that the membrane adsorber capsule can be used in bind and elute applications. As a result, polishing economy is improved and buffer consumption is reduced.



#### Comparison of Binding Capacity: Standard vs. Void-Volume-Optimized Capsule


As a result of the minimized void volume, the dynamic binding capacity of the Sartobind 8 mm capsules is increased by nearly 50%.

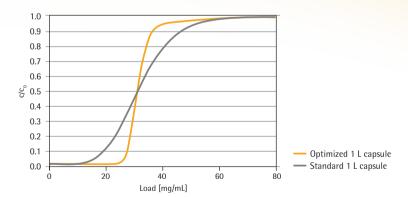

trypsimogen A

#### **Improved Resolution**

The efficient design also provides for sharp elution peaks and excellent resolution. The small peak | pool volume makes Sartobind membrane adsorbers applicable for bind and elute processes.

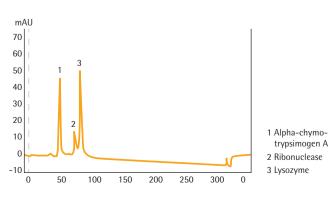

# For Virus | Protein Capture and Contaminant Removal

With macropores larger than 3  $\mu$ m, Sartobind membrane adsorbers are especially suited for the purification of large molecules, such as blood factors and virus particles.

They are also successfully used in flow-through polishing of recombinant proteins like monoclonal antibodies to remove host cell proteins, DNA, aggregates and other contaminants. These polishing steps are performed at high flow rates of up to 5 bed volumes per minute.



3 Determination Capacity

Optimizing Conditions in 96-Well Plates

Ligand Selection

#### Ligands Applications - Available Optimized Format

Membrane Type, Capacity	Single Use	Reuse	For Polishing	For Capture	4 mm	8 mm
Q anion exchanger 29 mg/mL with BSA	<b>✓</b>	<b>✓</b>	DNA, host cell proteins, endotoxins	Viruses, VLP, large proteins (e.g. FVIII)	<b>✓</b>	<b>✓</b>
S cation exchanger 25 mg/mL with lysozyme	<b>√</b>	<b>✓</b>	Host cell proteins, aggregates	Viruses, VLP, large proteins	<b>√</b>	<b>✓</b>
Phenyl hydrophobic inter- action 15 mg/mL with IgG	<b>√</b>	<b>√</b>	Aggregates, detergents, viruses	Viruses, VLP, large proteins	-	<b>√</b>

4



# of Binding with Sartobind pico or nano

## Accessories

Sartobind Jumbo 5 L on a trolley



Adapter for use with a 150 mL capsule on a liquid chromatography system






### Frequently Asked Questions

#### What is the direction of flow?

While the fluid enters at the top of the capsule, it flows from the external upstream to the internal downstream channel.

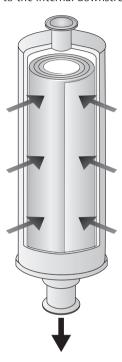

#### How is the membrane stacked?

The membranes are wheeled up, forming the membrane bed.

#### Can I reuse Sartobind capsules?

Yes, you can. The membrane type Q has been reused for 1,000 cycles while the HIC has already been reused successfully for 200 cycles.

# Can I test the integrity of the adsorbers before and after using them?

Yes. A diffusion test for this purpose is described in the manual.

#### Can I run capsules in series or parallel?

The capsules have been run successfully in both ways as well as in the combination of both.

## What type of device should I start with in the lab?

After selecting your ligand, trials can be run with Sartobind pico. At a 0.08 mL bed volume, it requires only a minimum sample quantity for binding studies. For screening binding conditions, 96-well plates can be used.

### How do I choose between using a column or a membrane adsorber?

Columns are recommended for bind and elute applications, where their capability of binding small molecules can be best employed. However, for rapid contaminant removal – especially when the process volume is large – and the capture of viruses or large proteins, Sartobind capsules are an excellent choice.

#### Can the capsules be autoclaved?

Yes, the ion exchange capsules can be autoclaved once at 121°C for 30 minutes.

# When was the first commercial process approved using Sartobind?

The monoclonal antibody Campath 1-H was approved in 2001. Therefore, Sartobind has been used for DNA removal in production ever since.

# Sartobind capsules are intended for single-use applications. Yet the literature states the capsules may be reused. How do I decide whether to reuse the capsules?

Typically, membrane adsorbers are used only once. Whether it makes sense for you to reuse them can be determined by cost-benefit calculations, which take the validation and cleaning costs of reuse into account.

### **Technical Data**

#### Membrane

Base material	Stabilized reinforced cellulose
Membrane thickness	275 μm
Pore size	> 3 µm
lon exchanger types	– Strong anion Q (quaternary ammonium) R-CH <sub>2</sub> -N <sup>+</sup> (CH <sub>3</sub> ) <sub>3</sub> – Strong cation S (sulfonic acid) R-CH <sub>2</sub> -SO $_3^-$
Hydrophobic interaction type	– Phenyl (-NH-C <sub>6</sub> H <sub>6</sub> )

#### **Capsule Materials**

Outer cage, inner core, end caps, capsule housings, nonwovens	Polypropylene
O-ring in vent valve*	Silicone

<sup>\*</sup>Only with 150 mL, 1.2 L and 5 L  $\,$ 

#### Operation

Depyrogenation	1 N NaOH for 30 minutes at 20°C
Autoclaving	121°C for 30 minutes for one cycle only (for Q and S)
Integrity testing	By the diffusion test method

#### **Dimensions and Connections**

Capsule size	nano 3 mL	75 mL   150 mL	1.2 L	5 L Jumbo
			ſ	
			100	1200
	ġ.	Ĭ.	Į.	
Bed height	8 mm	4 mm   8 mm	8 mm	8 mm
Membrane (cm²)	110	2700   5,500	44,000	182,000
Dimensions (height × diameter) mm	37×33	190×77	810×100	850×302
Connectors	Luer female	Sanitary 3/4"	Sanitary 1 1/2"	Sanitary 1 1/2"
		25 mm outer,	50.5 mm outer,	50.5 mm outer,
		14 mm inner diameter	36 mm inner diameter	36 mm inner diameter
Dimensions hose barb* (height × diameter) mm	-	203×77	-	-
Approximate weight	10 g	400 g	1.9 kg	16 kg, 20 kg wet, 23 kg filled

<sup>\*</sup> For 75 mL capsules only

Related products are Sartobind 96-well plates, MA units for laboratory use and the Sartobind capsules with a 4 mm bed height: Sartobind nano, 1 mL; mini, 7 mL; 5", 70 mL; 10", 180 mL; 20", 360 mL; 30", 540 mL; and mega, 1.62 L.

#### For more information:

- Sartobind brochure, Publication No.: SL-1513-e, Order No.: 85030-518-92
- Sartobind brochure for Polishing, Publication No.: SL-1521-e, Order No.: 85030-531-08

# **Ordering Information**

#### Sartobind Q Capsules

Order no.	Chemistry	Description adapter inlet and outlet	Quan- tity	Bed volume [mL]	Bed height [mm]	Nominal membrane area	Protein binding capacity** [g]	Recom- mended flow rate [L/min]	Approx. weight
Sartobind nano 96IEXQ42EUC11A	Q	Female Luer	4	3	8	110 cm <sup>2</sup>	0.088	0.015	10 g
Sartobind 75 mL 96IEXQ42D9M00A	Q	Hose barb	4	75	4	2,700 cm <sup>2</sup>	2.16	1.5	400 g
Sartobind 75 mL 96IEXQ42D9MFFA	Q	Sanitary flange	4	75	4	2,700 cm <sup>2</sup>	2.16	1.5	400 g
Sartobind 150 mL 96IEXQ42E9BFF	Q	Sanitary flange	1	150	8	5,500 cm <sup>2</sup>	4.4	0.75	400 g
Sartobind 1.2 L 96IEXQ42E3FSS	Q	Sanitary flange	1	1,200	8	4.4 m <sup>2</sup>	35	6	1.9 kg
Sartobind Jumbo 5 L 96IEXQ42E3ESS	Q	Sanitary flange	1	5,000	8	18.2 m <sup>2</sup>	45	25	16 kg
Sartobind S Capsules									
Order no.	Chemistry	Description adapter inlet and outlet	Quan- tity	Bed volume [mL]	Bed height [mm]	Nominal membrane area	Protein binding capacity** [g]	Recom- mended flow rate [L/min]	Approx. weight
Sartobind nano 96IEXS42EUC11A	S	Female Luer	4	3	8	110 cm <sup>2</sup>	0.077	0.015	10 g
Sartobind 75 mL 96IEXS42D9M00A	S	Hose barb	4	75	4	2,700 cm <sup>2</sup>	1.89	1.5	400 g
Sartobind 75 mL 96IEXS42D9MFFA	S	Sanitary flange	4	75	4	2,700 cm <sup>2</sup>	1.89	1.5	400 g
Sartobind 150 mL 96IEXS42E9BFF	S	Sanitary flange	1	150	8	5,500 cm <sup>2</sup>	3.9	0.75	400 g
Sartobind 1.2 L 96IEXS42E3FSS	S	Sanitary flange	1	1,200	8	4.4 m <sup>2</sup>	31	6	1.9 kg
Sartobind Jumbo 5 L 96IEXS42E3ESS	S	Sanitary flange	1	5,000	8	18.2 m <sup>2</sup>	127	25	16 kg
Sartobind Phenyl Caps	sules								
Order no.	Chemistry	Description adapter inlet and outlet	Quan- tity	Bed volume [mL]	Bed height [mm]	Nominal membrane area	Protein binding capacity** [g]	Recom- mended flow rate [L/min]	Approx. weight
Sartobind nano 96HICP42EUC11A	Phenyl	Female Luer	4	3	8	110 cm <sup>2</sup>	0.044	0.015	10 g
Sartobind 150 mL 96HICP42E9BFF	Phenyl	Sanitary flange	1	150	8	5,500 cm <sup>2</sup>	2.2	0.75	400 g

Phenyl

Phenyl

Sartobind 1.2 L

96HICP42E3FSS

96HICP42E3ESS

Sartobind Jumbo 5 L

1

1,200

5,000

8

8

 $4.4 \text{ m}^2$ 

18.2 m<sup>2</sup>

18

73

6

25

1.9 kg

16 kg

Sanitary

Sanitary

flange

flange

<sup>\*</sup> Frontal surface area = average cylinder surface

<sup>\*\*</sup> Typical dynamic binding capacity at 10% breakthrough was measured with BSA for Q | Iysozyme for S chemistry and IgG for HIC. 1 mL is the equivalent to 36 cm² membrane

#### Accessories

Order No.	Description	Quantity	
1ZA0GV0003	Adapter, Sanitary 25 mm, $\frac{1}{2}$ " - $\frac{3}{4}$ " to UNF 10 – 32	2	
1ZA0004	Adapter, male Luer to UNF 10-32 female, PEEK	1	
9ZGL0102	Trolley for Jumbo, 5 L, stainless steel	1	
7ZAL-V0013	Reducing adapter 1 ½" - ¾" / ¼"; 50.5 mm to 25 mm, sanitary	2	
7ZAL-V0010	Reducing adapter 2" - 1 1/2"; 64/50.5 mm to 25 mm, sanitary	2	

#### 96-Well Plates and Accessories

Order No.	Description	Quantity
99IEXQ42GCV	Sartobind Q 96-well plate, 2 units	2 (24 × 8-well strips)
99IEXQ42GCD	Sartobind Q 96-well plate, 10 units	10 (120 × 8-well strips)
99STPA42GCV	Sartobind STIC PA 96-well plate, 2 units	2 (24 × 8-well strips)
99STPA42GCD	Sartobind STIC PA 96-well plate, 10 units	10 (120 × 8-well strips)
99IEXS42GCV	Sartobind S 96-well plate, 2 units	2 (24 × 8-well strips)
99IEXS42GCD	Sartobind S 96-well plate, 10 units	10 (120 × 8-well strips)
99HICP42GCV	Sartobind Phenyl 96-well plate, 2 units	2 (24 × 8-well strips)
99HICP42GCD	Sartobind Phenyl 96-well plate, 10 units	10 (120 × 8-well strips)
VW96VAC01	Vac96 vacuum manifold	1
VW96VAA02	Vac96 liquid trap and reservoir	1
VW96VAA04	96-deep-well collection plate, 2 mL (square wells)	25
VW96VAA05	Replacement seal for Vac96 vacuum manifold	25
VW08VAC01	Vac8 vacuum manifold	1
VW08VAA02	Vac8 liquid trap and reservoir	1
VW08VAA03	8-well collection strips, 1.2 mL (round wells)	125
VW08VAA04	Replacement seal for Vac8 vacuum manifold	1
BBI-8582572	Swing-out rotor for deep-well plates (max. 80 mm high), incl. 2 buckets	1
16694-2-50-06	Microsart <sup>®</sup> mini.vac vacuum pump, 230 V∣50 Hz	1
16694-1-60-06	Microsart mini.vac vacuum pump, 115 V   60 Hz	1

#### **Sartobind pico Capsules**

Order No.	Description	Quantity
92IEXQ42DD11D	Sartobind Q pico 0.08 ml	10
92STPA42DD11D	Sartobind STIC PA pico 0.08 ml	10
92IEXS42DD11D	Sartobind S pico 0.08 ml	10
92HICP42DD11D	Sartobind Phenyl pico 0.08 ml	10

#### Sales and Service Contacts

For further contacts, visit www.sartorius-stedim.com

#### Europe

#### Germany

Sartorius Stedim Biotech GmbH August-Spindler-Strasse 11 37079 Goettingen

Phone +49.551.308.0 Fax +49.551.308.3289

Sartorius Stedim Systems GmbH Robert-Bosch-Strasse 5 – 7 34302 Guxhagen

Phone +49.5665.407 0 Fax +49.5665.407.2200

#### France

Sartorius Stedim FMT S.A.S. ZI des Paluds Avenue de Jouques - CS 91051 13781 Aubagne Cedex

Phone +33.442.845600 Fax +33.442.845619

Sartorius Stedim France SAS 71 des Paluds Avenue de Jouques - CS 71058 13781 Aubagne Cedex

Phone +33.442.845600 Fax +33.442.846545

#### Austria

Sartorius Stedim Austria GmbH Franzosengraben 12 1030 Vienna

Phone +43.1.7965763.18 Fax +43.1.796576344

Sartorius Stedim Belgium N.V. Leuvensesteenweg, 248/B 1800 Vilvoorde

Phone +32.2.756.06.80 Fax +32.2.756.06.81

#### Hungary

Sartorius Stedim Hungária Kft. Kagyló u. 5 2092 Budakeszi

Phone +36.23.457.227 Fax +36.23.457.147

#### Italy

Sartorius Stedim Italy S.p.A. Via dell'Antella, 76/A 50012 Antella-Bagno a Ripoli (FI)

Phone +39.055.63.40.41 Fax +39.055.63.40.526

#### Netherlands

Sartorius Stedim Netherlands B.V.

Phone +31.30.60.25.080 Fax +31 30 60 25 099

filtratie.nederland@sartorius-stedim.com

#### Poland

Sartorius Stedim Poland Sp. z o.o. ul Wrzesinska 70 62-025 Kostrzyn

Phone +48.61.647.38.40 Fax +48.61.879.25.04

#### Russian Federation

LLC "Sartorius ICR" Uralskaya str. 4, Lit. B 199155, Saint-Petersburg

Phone +7.812.327.5.327 Fax +7.812.327.5.323

Sartorius Stedim Spain, S.A.U. Avda. de la Industria, 32 Edificio PAYMA 28108 Alcobendas (Madrid)

Phone +34.902.110.935 Fax +34.91.358.96.23

#### Switzerland

Sartorius Stedim Switzerland AG Ringstrasse 24 a 8317 Tagelswangen

Phone +41.52.354.36.36 Fax +41.52.354.36.46

Sartorius Stedim UK Ltd. Longmead Business Centre Blenheim Road, Epsom Surrey KT19 9 QQ

Phone +44.1372.737159 Fax +44.1372.726171

#### America

Sartorius Stedim North America Inc. 5 Orville Drive, Suite 200 Bohemia NY 11716

Toll-Free +1.800.368.7178 Fax +1.631.254.4253

#### Argentina

Sartorius Argentina S.A. Int. A. Ávalos 4251 B1605ECS Munro **Buenos Aires** 

Phone +54.11.4721.0505 Fax +54.11.4762.2333

#### Brazil

Sartorius do Brasil Ltda Avenida Senador Verqueiro 2962 São Bernardo do Campo CEP 09600-000 - SP- Brasil

Phone +55.11.4362.8900 Fax + 55.11.4362.8901

#### Mexico

Sartorius de México S A de C V Circuito Circunvalación Poniente No. 149 Ciudad Satélite 53100, Estado de México México

Phone +52.5555.62.1102 Fax +52.5555.62.2942

#### Asia | Pacific

Sartorius Stedim Australia Pty. Ltd. Unit 5, 7-11 Rodeo Drive Dandenong South Vic 3175

Phone +61.3.8762.1800 Fax +61.3.8762.1828

Sartorius Stedim Biotech (Beijing) Co. Ltd. No. 33 Yu'an Road Airport Industrial Park Zone B Shunyi District, Beijing 101300

Phone +86.10.80426516 Fax +86.10.80426580

Sartorius Stedim Biotech (Beijing) Co. Ltd. Shanghai Branch Office 3rd Floor, North Wing, Tower 1 No. 4560 Jin Ke Road Pudong District, Shanghai 201210

Phone +86.21.68782300 Fax +86.21.68782332 | 68782882 Sartorius Stedim Biotech (Beijing) Co. Ltd. Guangzhou Representative Office Unit K, Building 23 Huihua Commerce & Trade Building No. 80 Xianlie Middle Road Guangzhou 510070

Phone +86.20.37618687 | 37618651 Fax +86.20.37619051

Sartorius Stedim India Pvt. Ltd. #69/2-69/3. NH 48. Jakkasandra Nelamangala Tg 562 123 Bangalore, India

Phone +91.80.4350.5250 Fax +91.80.4350.5253

#### Japan

Sartorius Stedim Japan K.K. 4th Fl., Daiwa Shinagawa North Bldg. 8-11, Kita-Shinagawa 1-chome Shinagawa-ku, Tokyo, 140-0001 Japan

Phone +81.3.4331.4300 Fax +81.3.4331.4301

#### Malaysia

Sartorius Stedim Malaysia Sdn. Bhd. Lot L3-E-3B, Enterprise 4 Technology Park Malaysia Bukit Jalil 57000 Kuala Lumpur, Malaysia

Phone +60.3.8996.0622 Fax +60.3.8996.0755

#### Singapore

Sartorius Stedim Singapore Pte. Ltd. 1 Science Park Road, The Capricorn, #05-08A Singapore Science Park II Singapore 117528

Phone +65.6872.3966 Fax +65.6778.2494

#### South Korea

Sartorius Korea Biotech Co., Ltd. 8th Floor, Solid Space B/D, PanGyoYeok-Ro 220, BunDang-Gu SeongNam-Si, GyeongGi-Do, 463-400

Phone +82.31.622.5700 Fax +82.31.622.5799



Chromatography as easy as filtration



