### **Instructions for Use**

# Vivaflow 50|50R|200



85030-519-25





## Contents

Vivaflow 50	3
Single Module	3
Linked Modules	4
Diafiltration   Desalting	5
Concentration	6
Recovery	6
System Components	7
Technical Specifications	7
Performance Characteristics	8
Ordering Information	9
Chemical Compatibility	10
Vivaflow 50R and Vivaflow 200	11
Operation of Single Module	11
Operation of Two Modules	12 – 13
Diafiltration   Desalting	14
Concentration	14
Recovery	14
Recovery Technical Specifications	14 15
· ·	
Technical Specifications	15
Technical Specifications Cleaning	15 16
Technical Specifications Cleaning Storage	15 16 16

#### Vivaflow 50

Designed for **single use** only. Anything else constitutes improper use.

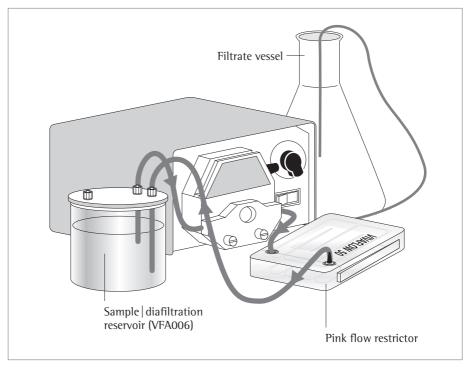


Diagram 1: Vivaflow 50 set up

#### Single module

- Set up the system as illustrated in Diagram 1. Note the positioning of the flow restrictor on the return line.
- Vivaflow membranes contain trace amounts of glycerine and sodium azide.
   To remove these chemicals and to check the security of the tube connections, it is recommended to rinse the module, and to test the system at full pressure before introducing the sample.
- 3. Place 500 ml deionised water in a suitable reservoir. The Vivaflow sample | diafiltration reservoir (prod. no. VFA006) is recommended.

- 4. Pump liquid through the system to purge any air pockets. The recirculation rate should be in the range 200–400 ml/min, and suitable flow should exit the filtrate line. If used, the pressure indicator (prod. no. VFA020), should read approximately 2.5 har.
- 5. Allow 400 ml to pass into the filtrate vessel. Check for any leakage at tubing connection points. Drain the system and empty or replace the filtrate vessel, (see recovery section). The system is now ready for use.

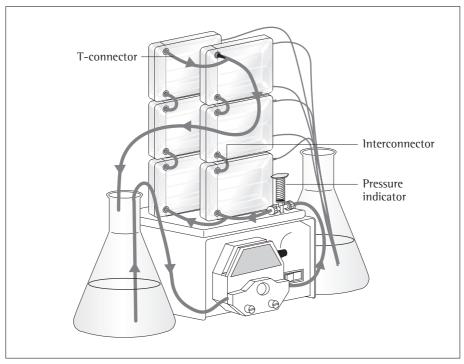


Diagram 2: Setting up multiple Vivaflow 50 cassettes

#### Linked modules

- Set up the system as illustrated in Diagram 2. Link the required number of modules by sliding the tongue and groove edges together. Note the positioning of the flow restrictor on the return line.
- Use series interconnectors, (prod. no. VFA031), to connect a single row of modules. Detailed ordering information for system components to be found of page 7, Diagram 4.
- 3. In addition, use T-connectors, (prod. no. VFA030), to connect two rows of modules in parallel.
- 4. The flow path of the system can be configured to suit the membrane | sample combination. Most solutions are better suited by modules connected in parallel. Very low viscosity solutions favour modules connected only in series. Rinse the system as detailed under single module operation.

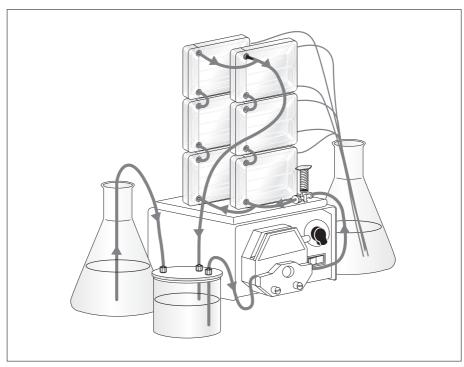


Diagram 3: Buffer exchange | diafiltration with Vivaflow 50

#### Diafiltration Desalting

- Set up the system as illustrated in Diagram 3 for linked modules. If only one module is required, connect it in place of the linked modules.
- 2. Rinse the system as detailed previously.
- 3. Either place the solution to be purified in the 500 ml sample | diafiltration reservoir (prod. no. VFA006), or concentrate a larger volume to 500 ml or less so that it is fully contained in the 500 ml reservoir. Ensure the reservoir lid is firmly closed.
- Fill a large feed reservoir with exchange solvent and pump sample through the system as for concentration.

5. As the volume in the 500 ml reservoir decreases, the vacuum created draws exchange solvent through the feed line from the larger reservoir. Over 99% solvent exchange can be accomplished with an exchange volume approximately 5 times the volume of the sample.

#### Concentration

- 1. Fill the feed reservoir with sample solution. When initial volumes larger than 500 ml are required, place 500 ml in the sample | diafiltration reservoir, the remaining volume in another suitable container and connect the vessels as detailed in the diafiltration section. Alternatively, use a larger container for the entire sample volume and immerse the feed and return lines directly into the liquid.
- 2. Pump liquid through the system. The recirculation rate should be in the range 200–400 ml/min, and suitable flow should exit the filtrate line. If used, the pressure indicator, (prod. no. VFA020), should read approximately 2.5 bar.
- 3. Concentrate the sample to the desired volume.
- ▲ Warning: Do not run the same section of tubing through the pump head for longer than six hours, over use of tubing will result in significant pressure drop and ultimately, failure.
- 4. When the desired volume has been reached, reduce the recirculation rate to 20–40 ml/min and recirculate the concentrated sample for 1–2 minutes to maximise recovery.

#### Recovery

- 1. Disconnect the feed line from the lid of the 500 ml reservoir or when using a different container, remove the feed line from the sample.
- Pump residual system volume back into the reservoir | container. (When parallel modules are used with viscous solutions, ensure that all modules are empty by pinching the tubing between each of the stacks of modules in turn).
- For a more complete sample recovery rinse approximately 5–10 ml per module of water or sample buffer through the system, and recover as before.

#### **System Components**

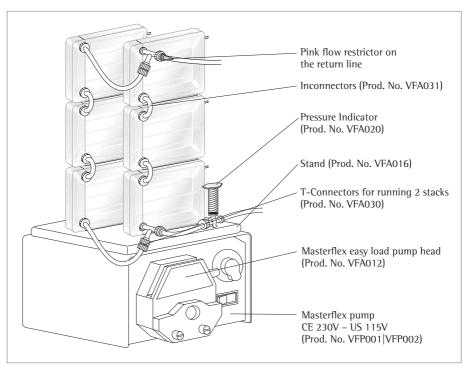


Diagram 4: System componets for setting up multiple Vivaflow 50 units

**Table 1: Technical Specifications** 

Dir	ne	nsı	n	ns

Overall L H W	107   84   25 mm
Channel W   H	15 mm   0.3 mm
Active membrane area	50 cm <sup>2</sup>
Hold up volume (module)	1.5 ml
Min recirculation volume	< 10 ml
Non recoverable hold-up	< 0.5 ml

#### **Operating Conditions**

Pump flow	200-400 ml/min
Maximum pressure	3 bar (45 psi)
Maximum temperature	60°C

#### **Materials of Construction**

Main housing	Polycarbonate
Flow channel	TPX (PMP)
Membrane support	TPX (PMP)
Seals and 0 rings	Silicone
Pressure indicator	Polypropylene, SS Spring,
Flow restrictor	Polypropylene
Fittings	Nylon
Tubing	PVC (medical grade)

**Table 2: Performance characteristics** 

		Time to concentrate up to 20x (min.) at 3 bar inlet pressure, 20°C			
	Single Device	Three Devices	Solute Recovery %		
	250 ml Start Vol. min.	1 L Start Vol. min.	Direct	10 ml rinse	
BSA 1.0 mg/ml (66,000 MW)	)				
5,000 MWCO PES	34	49	96%	> 99%	
10,000 MWCO PES	22	32	94%	> 99%	
30,000 MWCO PES	22	32	92%	99%	
50,000 MWCO PES	20	29	92%	98%	
γ Globulins 1.0 mg/ml					
100,000 MWCO PES	43	62	92%	98%	
100,000 MWCO RC	40	58	92%	98%	
100,000 MWCO Hydrosart	40	58	92%	98%	
Yeast 1.0 mg/ml (S.Cerevisia	e)				
0.2 um PES	33	47	92%	98%	

## **Ordering Information**

Vivaflow 50 include filtrate tube, size 16 peristaltic tubing, flow restrictor and fittings and $1 \times Series$ Interconnector	Pack Size	Prod. No.
3,000 MWC0 (PES)	2	VF05P9
5,000 MWC0 (PES)	2	VF05P1
10,000 MWCO (PES)	2	VF05P0
30,000 MWC0 (PES)	2	VF05P2
50,000 MWC0 (PES)	2	VF05P3
100,000 MWC0 (PES)	2	VF05P4
0.2 μm (PES)	2	VF05P7
100,000 MWCO (RC)	2	VF05C4
Vivaflow 50 complete system comprises		
Pump (240 V), easy load pump head (size 16), tubing, 500 ml sample   diafiltration reservoir, module stand, pressure indicator, T connectors, series interconnectors	1	VFS502
Pump (115 V), easy load pump head (size 16), tubing, 500 ml sample   diafiltration reservoir, module stand, pressure indicator, T connectors, series interconnectors  Vivaflow 50 PVC tubing and fittings	1	VFS504
Size 16 PVC pump tubing (3 metres, 3.2×1.6 mm)		VFA004
Flow restrictor set $(2 \times 0.4, 0.6, 0.8 \text{ mm})$		VFA009
T connectors for running 2 stacks (2 pieces)		VFA030
Series interconnectors (6 pieces)		VFA031
Female luer fittings (10 pieces)		VFA032
VF50 tubing kit (2×1 m size 16 PVC tubing with inlets fittings, 2×50 cm size 16 PVC tubing with 0.6 mm flow restrictors, 1× series inlet connections)		VFA034
Flow restrictor 0.6 mm (6 pieces)		VFA035
Vivaflow 50 accessories		
Masterflex economy drive variable speed peristaltic pump (230 V)		VFP001
Masterflex economy drive variable speed peristaltic pump (115 V)		VFP002
500 ml sample and or diafiltration reservoir		VFA006
Masterflex easy load pump head – size 15		VFA013
Masterflex easy load pump head – size 16		VFA012
Vivaflow 50 stand		VFA016
Pressure indicator (1–3 bar)		VFA020

Table 3: Chemical Compatibility for the Vivaflow 50

Solution	PES	RC
Acetic acid (25%)	OK	OK
Acetone	NO	NO
Ammonium hydroxide (5%)	OK	OK
Ammonium sulfate sat.	OK	OK
DMEM	OK	OK
Ethanol (70%)	OK	OK
Ethyl acetate	NO	NO
Formaldehyde (30%)	OK	OK
Formic acid (5%)	OK	OK
Guanidine HCl (6 M)	OK	OK
Hydrocarbons, Aromatic	NO	?
Hydrocarbons, Chlorinated	NO	NO
Hydrochloric acid (1 M)	OK	NO
Isopropanol	NO	NO
Lactic acid (5%)	OK	OK
Mercaptoethanol (10 mM)	OK	OK
Methanol (60%)	OK	OK
n-Butanol (70%)	OK	OK

Solution	PES	RC
Peracetic acid (0.2%)	OK	OK
Phenol (1%)	OK	?
Phosphate Buffer (1 M)	OK	OK
Pyridine	NO	NO
RPMI-1640	OK	OK
Sodium azide	OK	OK
Sodium deoxycholate (5%)	OK	OK
Sodium hydroxide (1 M)	NO	NO
Sodium hydroxide (0.1 M)	NO	NO
Sodium hypochlorite (0.02%)	?	?
Sodium nitrate	NO	NO
Sulfamic acid (5%)	OK	NO
Surfactants (0.1%)	OK	OK
Toluene	NO	NO
Trichloroacetic acid (10%)	NO	?
Trifluoroacetic acid (10%)	OK	NO
Urea (8 M)	OK	OK

OK = Acceptable ? = Questionable NO = Not recommended

#### Vivaflow 50R and Vivaflow 200

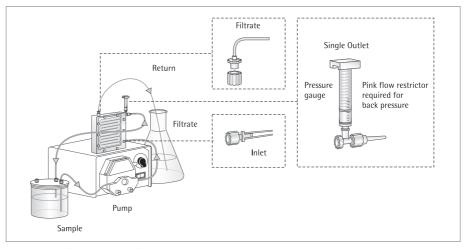


Diagram 5: Vivaflow 50R and Vivaflow 200 set up

## Operation Single Vivaflow 50R or Vivaflow 200 Modules

#### Assembly and rinsing single module

- 1. Set up the system as illustrated above in Diagram 5.. Note the positioning of the flow restrictor on the return line. A set of flow restrictors of different gauges (Product No. VFA009), is available for use with solutions of unusually high or low viscosity.
- Vivaflow membranes contain trace amounts of glycerine and sodium azide. To remove these chemicals and to check the security of the tube connections, it is recommended to rinse the module, and to test the system at full pressure before introducing the sample.
- 3. Place 500 ml deionised water in a suitable reservoir.

- 4. Pump liquid through the system to purge any air pockets. The recirculation rate should be 200–400 ml/min, and suitable flow should exit the filtrate line. If used, the pressure indicator should read approximately 2.5 bar.
- Allow 400 ml to pass into the filtrate vessel. Check for any leakage at tubing connection points.
- Drain the system and empty or replace the filtrate vessel (see Recovery section). The system is now ready for use.

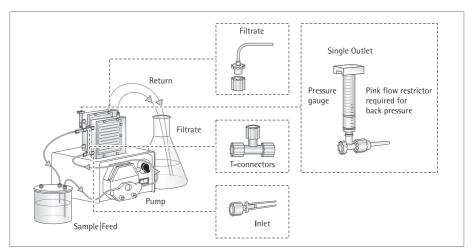


Diagram 6: Setting up two Vivaflow 50R units

## Operating two Vivaflow 50R or two Vivaflow 200 Modules

#### Assembly and rinsing two modules

1. Set up the Vivaflow 50R system as illustrated in Diagram 6 and the Vivaflow 200 System as illustrated in Diagram 7 system as illustrated above.

#### 2. Vivaflow 50R:

Use two T-connectors (VFA030) to connect to the inlets of both Vivaflow 50R modules.

#### Vivaflow 200:

Use the Y-connector, (VFA005) to connect to the inlets of both Vivaflow 200 modules.

The size 15 tubing connected to the Y-connector is the inlet tubing for the pump. Use a size 15 pump head.

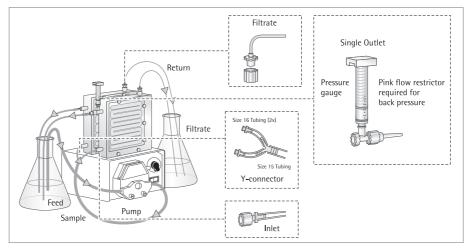


Diagram 7: Setting up two Vivaflow 200 units

- 3. Note the positioning of a flow restrictor on the return line from each Vivaflow 200 module. Vivaflow membranes contain trace amounts of glycerine and sodium azide. To remove these chemicals and to check the security of the tube connections, it is recommended to rinse the module, and to test the system at full pressure before introducing the sample.
- 4. Place 1 litre of deionised water in a suitable reservoir.
- 5. Pump liquid through the system to purge any air pockets. The recirculation rate should be 200-400 mL/min (500-900 mL/min for two Vivaflow 200 cassettes), and suitable flow should exit the filtrate line. If used, the pressure indicator should read approximately 2.5 bar.

- 6. Allow 800 ml to pass into the filtrate vessel. Check for any leakage at tubing connection points.
- 7. Drain the system and empty or replace the filtrate vessel (see Recovery section). The system is now ready for use.

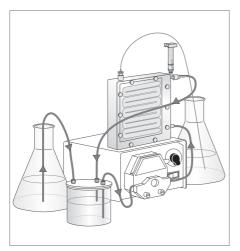


Diagramm 8: Buffer exchange | Desalting with Vivaflow 50R and Vivafow 200

#### Diafiltration Desalting

- 1. Set up the system as illustrated in Diagram 8.
- 2. Rinse the system as detailed previously.
- 3. Place the solution to be purified in a suitable diafiltration reservoir. For volumes up to 500 ml the Vivaflow 500 ml sample diafiltration reservoir (VFA006), is recommended. Ensure the lid is firmly closed.
- Fill the large feed reservoir with exchange solvent and pump sample through the system as for concentration.
- 5. As liquid passes through the membrane, the vacuum created in the diafiltration reservoir draws in exchange solvent through the feed line from the large reservoir. Over 99% solvent exchange can be accomplished with an exchange volume of approximately 5 times the sample volume.

#### Concentration

- 1. Fill the feed reservoir with sample solution.
- Pump liquid through the system.
   The recirculation rate should be 200–400 ml/min, (500–900 ml/min for two modules), and suitable flow should exit the filtrate line. If used, the pressure indicator should read approximately 2.5 bar.
- ▲ Warning: Do not run the same section of tubing through the pump-head for longer than 6 hours, overuse of tubing will result in significant pressure drop and ultimately, failure.
- 3. Concentrate the sample.
- When nearing the desired volume, reduce the recirculation rate to 20–40 ml/min and recirculate the concentrated sample for 1–2 minutes to increase sample recovery.

#### Recovery

- 1. Remove the feed line from the sample.
- 2. Pump residual system volume back into the reservoir container.
- 3. For a more complete recovery, rinse 25–50 ml of water or sample buffer through the system, and recover as before.

Table 4: Technical specifications Vivaflow 50R and Vivaflow 200

Dimensions	Vivaflow 50 R	Vivaflow 200	
Overall L H W	100   100   24 mm	126   138   38 mm	
Channel W   H	7.5   0.4 mm	10 mm   0.4 mm	
Active membrane area	50 cm <sup>2</sup>	200 cm <sup>2</sup>	
Hold up volume (module)	1.7 ml	5.3 ml	
Min. recirculation volume	10 ml	< 20 ml	
Non recoverable hold-up	< 0.5 ml	< 1 ml	
Materials of construction	Vivaflow 50R and Vivaflow	200	
Main housing	Acrylic		
Flow channel	Acrylic		
Membrane support	Polypropylene		
Seals and O rings	Silicone		
Pressure indicator	Polypropylene, SS spring		
Flow restrictor	Polypropylene		
Fittings	Nylon		
Tubing	PVC (medical grade)		
Operating conditions	Vivaflow 50R and Vivaflow	200	
Pump flow	200-400 ml/min		
Maximum pressure	4 bar (60 psi)		
	Pressure drop across inlet	outlet 0.5 bar (7 psi)	
Maximum temperature	60°C		

# Cleaning and storage information only for Vivaflow 50R and Vivaflow 200

#### Cleaning

Vivaflow 50R and 200 modules may be used several times if cleaning and storage instructions are followed.

- Flush the system with 200 ml of deionised water with the filtrate going to waste.
- 2. Place the feed, return and filtrate lines in a suitable container for the cleaning solution.
- 3. Prepare cleaning solutions suitable for the membrane.
- Polyethersulfone membranes:
   250 ml of 0.5 mM NaOCl in 0.5 M NaOH
- II. Hydrosart membranes: 250 ml of 0.5 M NaOH
- 4. Recirculate at 50–100 ml/min for 30–40 minutes.
- Drain the system and recirculate 250 ml of deionised water through the system for 5–10 minutes.
- 6. Drain and discard rinse solution and flush with a further 500 ml of deionised water with the filtrate going to waste. The system is now ready for further use.

#### Storage

To store Vivaflow 50R and Vivaflow 200 after cleaning, fill module with deionised water and 10% ethanol. Seal inlet, outlet and filtrate ports and refrigerate at approximately 4°C.

**Table 5: Performance characteristics Vivaflow 50R** 

Time to concentrate up to 20x (min) at 3.0 bar inlet | 2.5 bar outlet pressure, 20°C

	start volume	Average flux	Recovery %	0
	250 ml ml/min	Direct	25 ml rinse	
Lysozyme 0.25 mg/ml (14,000	MW)			
5,000 MWCO Hydrosart	70	3.4	96%	98%
10,000 MWCO Hydrosart	23	10.3	94%	96%
BSA 1.0 mg/ml (66,000 MW)				
10,000 MWCO Hydrosart	24	9.9	98%	> 99%
30,000 MWCO Hydrosart	15	15.8	97%	> 99%
γ Globulins 1.0 mg/ml (150,000	0 MW)			
100,000 MWCO Hydrosart	46	5.2	97%	> 99%
Start volume 1 litre (one Vivaf	low 50R at 3 bar)	10,000 MWC0	) Hydrosart	
BSA 1.0 mg/ml	95	10.0	98%	> 99%
Start volume 1 litre (two Vivaf	low 50R in paralle	el at 3 bar) 10,	000 MWCO	Hydrosart
BSA 1.0 mg/ml	48	19.8	98%	> 99%

**Table 6: Performance characteristics Vivaflow 200** 

Time to concentrate up to 20x (min.) at 3 bar inlet pressure, 20°C

	start volume	Average flux	Recovery	%
	1 litre	ml/min	Direct	25 ml rinse
Lysozyme 0.25 mg/ml (14,000	MW)			
2,000 MWCO Hydrosart	160	6	97%	> 99%
3,000 MWCO PES	180	5	97%	> 99%
BSA 1.0 mg/ml (66,000 MW)				
5,000 MWCO PES	29	33	98%	> 99%
5,000 MWCO Hydrosart	70	14	98%	> 99%
10,000 MWCO PES	23	41	96%	> 99%
10,000 MWCO Hydrosart	35	27	98%	> 99%
30,000 MWCO PES	25	38	96%	99%
30,000 MWCO Hydrosart	20	48	96%	> 99%
50,000 MWCO PES	22	43	96%	98%
$\gamma$ Globulins 1.0 mg/ml (averag	e 160,000 MW)			
100,000 MWCO PES	54	18	96%	99%
Yeast 1.0 mg/ml (S. Cerevisiae	e)			
0.2 μm PES	11	86	92%	98%
Dilute solute concentration, s	tart volume 1 litr	e at 3 bar, 10,0	00 MWCO	PES
BSA 0.001 mg/ml	18	52	90%	98%
BSA 0.01 mg/ml	20	47	92%	98%
BSA 0.1 mg/ml	21	45	94%	99%
Start volume 5 litres (two Viv	aflow 200 in para	allel at 3 bar) 10	0,000 MW	CO PES
BSA 1.0 mg/ml (66,000 MW)	67	70	97%	> 99%

## **Ordering Information**

Vivaflow 50R modules include pressure indicator, flow restrictor and size 16 pvc peristaltic tubing and fittings	Oty per box	Prod. no.
5,000 MWCO Hydrosart	1	VF05H1
10,000 MWCO Hydrosart	1	VF05H0
30,000 MWCO Hydrosart	1	VF05H2
100,000 MWCO Hydrosart	1	VF05H4
Vivaflow 200 modules include pressure indicator, flow restrictor and size 16 pvc peristaltic tubing and fittings	Oty per box	Prod. no.
3,000 MWCO PES	1	VF20P9
5,000 MWCO PES	1	VF20P1
10,000 MWCO PES	1	VF20P0
30,000 MWCO PES	1	VF20P2
50,000 MWCO PES	1	VF20P3
100,000 MWCO PES	1	VF20P4
0.2 μm PES	1	VF20P7
2,000 MWCO Hydrosart	1	VF20H9
10,000 MWCO Hydrosart	1	VF20H0
30,000 MWCO Hydrosart	1	VF20H2
100,000 MWCO Hydrosart	1	VF20H4
Vivaflow 50R   200 complete system comprises		
Pump (240 V), Easy load pump head (size 16), tubing, 500 ml sample   diafiltration reservoir	1	VFS202
Pump (115 V), Easy load pump head (size 16), tubing, 500 ml sample   diafiltration reservoir	1	VFS204

### Vivaflow 50R | 200 accessories

Masterflex economy drive variable speed peristaltic pump (230V)	VFP001
Masterflex economy drive variable speed peristaltic pump (115V)	VFP002
500 ml sample and or diafiltration reservoir	VFA006
Masterflex easy load pump head – size 16	VFA012
Masterflex easy load pump head – size 15 (only for operating two Vivaflow 200)	VFA013
Size 15 pvc pump tubing and Luer fittings (3 m, $4.8 \times 2.6$ mm)	VFA003
Size 16 pvc pump tubing and Luer fittings (3 m, $3.2 \times 1.6$ mm)	VFA004
Flow restrictor set $(2 \times 0.4, 0.6, 0.8 \text{ mm})$	VFA009
Female luer fittings size 16 (10 pieces)	VFA032
Flow restrictors 0.6 mm (6 pieces)	VFA035
Female luer fittings size 15 (10 pieces)	VFA036
Flow restrictor set (2 + 0.4, 0.6, 0.8 mm)	VFA009
Female luer fittings size 16 (10 pieces)	VFA032
Flow restrictors 0.6 mm (6 pieces)	VFA035
Female luer fittings size 15 (10 pieces)	VFA036
Vivaflow 200 tubing and fittings for operation two modules	
Y connector (size 15 to $2 \times$ size 16, Luer fittings)	VFA005
Masterflex easy load pump head – size 15	VFA013
Vivaflow 50R tubing a fittings for operation of two modules	
T-Connector	VFA030

Table 7: Chemical Compatibility for Vivaflow 50R and Vivaflow 200

Solution	PES	HY
Acetic acid (25%)	OK	OK
Acetone	NO	NO
Ammonium hydroxide (5%)	OK	OK
Ammonium sulfate sat.	OK	?
DMEM	OK	OK
Ethanol (70%)	OK	OK
Ethyl acetate	NO	NO
Formaldehyde (30%)	OK	OK
Formic acid (5%)	OK	OK
Guanidine HCl (6 M)	OK	OK
Hydrocarbons, Aromatic	NO	NO
Hydrocarbons, Chlorinated	NO	NO
Hydrochloric acid (1 M)	OK	OK
Isopropanol	NO	NO
Lactic acid (5%)	OK	OK
Mercaptoethanol (10 mM)	OK	OK
Methanol (60%)	OK	OK
n-Butanol (70%)	OK	OK

Solution	PES	HY
Peracetic acid (0.2%)	OK	NO
Phenol (1%)	OK	NO
Phosphate Buffer (1 M)	OK	OK
Pyridine	NO	NO
RPMI-1640	OK	OK
Sodium azide	OK	OK
Sodium deoxycholate (5%)	OK	?
Sodium hydroxide (1 M)	OK	OK
Sodium hydroxide (0.1 M)	OK	OK
Sodium hypochlorite (0.02%)	OK	NO
Sodium nitrate	NO	OK
Sulfamic acid (5%)	OK	?
Surfactants (0.1%)	OK	OK
Toluene	NO	NO
Trichloroacetic acid (10%)	NO	OK
Trifluoroacetic acid (10%)	OK	OK
Urea (8 M)	OK	OK

OK = Acceptable ? = Questionable

NO = Not recommended

Sartorius Stedim Lab Ltd. Sperry Way, Stonehouse Park GL10 3UT Stonehouse, Gloucestershire, UK

Phone: +44 1453 821972 www.sartorius.com

The information and figures contained in these instructions correspond to the version date specified below.

Sartorius reserves the right to make changes to the technology, features, specifications and design of the equipment without notice.

Masculine or feminine forms are used to facilitate legibility in these instructions and always simultaneously denote the other gender as well.

#### Copyright notice:

This instruction manual, including all of its components, is protected by copyright.

Any use beyond the limits of the copyright law is not permitted without our approval.

This applies in particular to reprinting, translation and editing irrespective of the type of media used.

Last updated: 08 | 2021

© 2021 Sartorius Stedim Lab Ltd. Sperry Way, Stonehouse Park GL10 3UT Stonehouse, Gloucestershire, UK

AM | Publication No.: SLU6097-e210810